



**STANDARD OPERATING PROCEDURE**

<b>Department:</b> Quality Control	<b>SOP No.:</b>
<b>Title:</b> Operation of Flask shaker	<b>Effective Date:</b>
<b>Supersedes:</b> Nil	<b>Review Date:</b>
<b>Issue Date:</b>	<b>Page No.:</b>

**1.0 OBJECTIVE:**

To lay down a procedure for operation of Flask shaker wrist action.

**2.0 SCOPE:**

This SOP is applicable for operation of Flask shaker wrist action in the Quality control department.

**3.0 RESPONSIBILITY:**

Officer, Executive – Quality Control.

Head– Quality Control.

**4.0 PROCEDURE:**

**4.1 Operation:**

4.1.1 Ensure that instrument is clean and free from dust.

4.1.2 Switch “ON” the mains.

4.1.3 Tighten the flask/beaker with the clamp and ensure that the flask/beaker is fitted with clamp.

4.1.4 Switch ON red power button of the instrument.

4.1.5 Adjust the shaking with black knob as per requirement. After completion of shaking switch OFF the instrument by the red power button.

**4.2 Calibration:**

4.2.1 Calibration of the instrument shall be done by engineering department.

**4.3 Precautions:**

4.3.1 Do not overload the clamp with too many samples.

4.3.2 Cleaning should be done with lint free duster frequently.

**5.0 ANNEXURE (S) :**

Nil



**PHARMA DEVILS**  
QUALITY CONTROL DEPARTMENT

**STANDARD OPERATING PROCEDURE**

<b>Department:</b> Quality Control	<b>SOP No.:</b>
<b>Title:</b> Operation of Flask shaker	<b>Effective Date:</b>
<b>Supersedes:</b> Nil	<b>Review Date:</b>
<b>Issue Date:</b>	<b>Page No.:</b>

**6.0 REFERENCE (S):**

SOP: Preparation, Approval, Distribution control, revision and destruction of Standard operating Procedure (SOP).

**7.0 ABBREVIATION (S)/DEFINITION (S):**

Nil

**REVISION CARD**

S.No.	REVISION No.	REVISION DATE	DETAILS OF REVISION	REASON (S) FOR REVISION	REFERENCE CHANGE CONTROL No.
1	00	---	---	New SOP	-