



**STANDARD OPERATING PROCEDURE**

<b>Department:</b> Quality Control	<b>SOP No.:</b>
<b>Title:</b> Operation of Heating Mantle	<b>Effective Date:</b>
<b>Supersedes:</b> Nil	<b>Review Date:</b>
<b>Issue Date:</b>	<b>Page No.:</b>

**1.0 OBJECTIVE:**

To lay down a procedure for Operation of Heating mental.

**2.0 SCOPE:**

This SOP is applicable for the use of Heating mental Mantle in the Quality Control Department.

**3.0 RESPONSIBILITY:**

Executive, Officer, Head of Department – Quality Control.

**4.0 PROCEDURE:**

**4.1 Operation:**

- 4.1.1 Ensure that instrument is clean and free from dust.
- 4.1.2 Switch ON the mains.
- 4.1.3 Adjust the thermostat to the required temperature by moving it in a clockwise direction.
- 4.1.4 Keep the vessels containing solution on the heating mantle.
- 4.1.5 After completion of work turn the knob to zero and switch OFF the main switch.

**4.2 Precaution:**

- 4.2.1 Do not handle flammable material near the instrument.
- 4.2.2 Clean the instrument properly after use.

**5.0 ANNEXURE (S):**

Nil

**6.0 REFERENCE (S):**

SOP: Preparation, Approval, Distribution control, revision and destruction of Standard Operating Procedure (SOP).

**7.0 ABBREVIATION (S)/DEFINITION (S):**

Nil



**PHARMA DEVILS**  
QUALITY CONTROL DEPARTMENT

**STANDARD OPERATING PROCEDURE**

<b>Department:</b> Quality Control	<b>SOP No.:</b>
<b>Title:</b> Operation of Heating Mantle	<b>Effective Date:</b>
<b>Supersedes:</b> Nil	<b>Review Date:</b>
<b>Issue Date:</b>	<b>Page No.:</b>

**REVISION CARD**

S. No.	REVISION No.	REVISION DATE	DETAILS OF REVISION	REASON (S) FOR REVISION	REFERENCE CHANGE CONTROL No.
1.	00	-	-	New Sop	-