

# PHARMA DEVILS

ENGINEERING DEPARTMENT

Title: Preventive Maintenance of Conta Blender

SOP No.:	Revision No.:	00
Effective Date:	Supersedes No.	Nil
Review Date:	Page No.	1 of 3

#### 1.0 OBJECTIVE

1.1 To describe a procedure for the preventive maintenance of Conta Blender.

#### 2.0 SCOPE

2.1 This procedure applies to the Engineering department.

#### 3.0 RESPONSIBILITY

3.1 Engineering

#### 4.0 ACCOUNTABILITY

4.1 Plant Head

### 5.0 REFERENCE (S)

5.1 In-house.

#### 6.0 PROCEDURE

- 6.1 Procedure for mechanical work:
  - 6.1.1 Switch off the machine and put "under maintenance" label.
  - 6.1.2 Clean the hydraulic tank from outside and tighten the connecting ferrules.
  - 6.1.3 Check the level of the Hydraulic oil in the tank if it is less than top up by servo 68.
  - 6.1.4 Clean the side plates and cover plates.
  - 6.1.5 Check the rotary movement of the machine.
  - 6.1.6 Grease the track from which the rollers moving.
- 6.2 Procedure for electrical work:
  - 6.2.1 Clean the panel board by hand blower.
  - 6.2.2 Clean the contactor and tighten the wire connection.
  - 6.2.3 Check the NO/NC element of UP/DOWN movement.
  - 6.2.4 Check the operation of limit switch.
  - 6.2.5 Check the alignment & direction of the motor.
  - 6.2.6 Check the connections of PLC.
- 6.3 Procedure for trial run:
  - 6.3.1 Switch on the power supply.
  - 6.3.2 Check the operation of the lifter referring the SOP.
  - 6.3.6 Check & record the observations as per Annexure-1.
- 6.4 Procedure for handover of the machines:
  - 6.4.1 Remove "under maintenance" label & affix the preventive maintenance status tag and inform the concerned department.
- Frequency of preventive maintenance is monthly.

#### 7.0 HISTORY

7.1 Details are given below.



# PHARMA DEVILS

ENGINEERING DEPARTMENT

Title: Preventive Maintenance of Conta Blender

SOP No.:	Revision No.:	00
<b>Effective Date:</b>	Supersedes No.	Nil
Review Date:	Page No.	2 of 3

SOP No.	REASON FOR CHANGE	EFFECTIVE DATE

# **8.0 ABBREVIATIONS**: The abbreviations used in the SOP are as follows:

8.1 SOP - Standard Operating Procedure

8.2 No. - Number

8.3 QA - Quality Assurance





ENGINEERING DEPARTMENT

Title: Preventive Maintenance	of Conta Blender		
SOP No.:		<b>Revision No.:</b>	00
<b>Effective Date:</b>		Supersedes No.	Nil
<b>Review Date:</b>		Page No.	3 of 3

# ANNEXURE-I

	Cleaning of the machine.	
_	Cicaming of the machine.	
2	Checking of operative units.	
3	Checking electrical connections.	
	Checking of wire rope & pulley alignment	
	Oil level in power pack.	