

# PHARMA DEVILS

ENGINEERING DEPARTMENT

<b>Title:</b> Preventive Maintenance of Infra Red Dryer
---

SOP No.:	Re	evision No.:	00
Effective Date:	Su	persedes No.	Nil
Review Date:	Pa	ge No.	1 of 3

#### 1.0 OBJECTIVE

1.1 To describe a procedure for the preventive maintenance of infra red dryer.

#### 2.0 SCOPE

**2.1** This procedure applies to the Engineering department.

#### 3.0 RESPONSIBILITY

3.1 Engineering

#### 4.0 ACCOUNTABILITY

4.1 Plant Head

### 5.0 REFERENCE (S)

5.1 In-house.

#### 6.0 PROCEDURE

- 6.1 Procedure for mechanical work:
  - 6.1.1 Switch off the power supply to the machine and put "under maintenance" label.
  - 6.1.2 Check for the wear and tear of PU coated castor wheels. If required, replace.
  - 6.1.3 Check for the smooth UP / DOWN movement of the umbrella; grease the guide rails/rollers assembly; if required.
  - 6.1.4 Clean the filter as per respective SOP.
- 6.2 Procedure for electrical work:
  - 6.2.1 Tighten the loose wire and connection of plug.
  - 6.2.2 Clean the operation of the controller.
  - 6.2.3 Clean the surface of heater (IR) by lint free cloth.
  - 6.2.4 Check heater bank for loose electrical connections.
- 6.3 Procedure for trial run:
  - 6.3.1 Switch on the power supply.
  - 6.3.2 Take the trial of machine by running the machine as per SOP of operation.
  - 6.3.3 Check & record the observations as per Format.
- 6.4 Procedure for handover of the machines:
  - 6.4.1 Remove under maintenance label & affix preventive maintenance status tag and inform the concerned department.
- 6.5 Frequency of preventive maintenance is once in two months.

#### 7.0 HISTORY

7.1 Details are given below.

SOP No.	REASON FOR CHANGE	EFFECTIVE DATE





ENGINEERING DEPARTMENT

Title: Preventive Maintenance of Infra Red Dryer

SOP No.:	Revis	sion No.:	00
Effective Date:	Super	rsedes No.	Nil
Review Date:	Page :	No.	2 of 3

**8.0 ABBREVIATIONS**: The abbreviations used in the SOP are as follows:

8.1 SOP - Standard Operating Procedure

8.2 No. - Number

8.3 QA - Quality Assurance





ENGINEERING DEPARTMENT

<b>Title:</b> Preventive Maintenance	of Infra Red Dryer		
SOP No.:		<b>Revision No.:</b>	00
<b>Effective Date:</b>		Supersedes No.	Nil
<b>Review Date:</b>		Page No.	3 of 3

## ANNEXURE-I

S.No.	Particulars	Done/Not Done	Remarks
1	Cleaning of machine.		
2	Check for the smooth functioning of the machine.		
3	Check heater bank for loose electrical connection.		
4	Check blower for loose electrical connections.		
5	Clean the filter.		
6	Cleaning of PU coated wheels and their brackets.		
7	Check all the fasteners.		
8	Check all the electrical connections.		