



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

**PERFORMANCE QUALIFICATION
REPORT
FOR
VIAL WASHING MACHINE**

EQUIPMENT ID. No.	
LOCATION	Washing & Sterilizing Room
DATE OF QUALIFICATION	
SUPERSEDES PROTOCOL No.	NIL



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

REPORT CONTENTS

S.No.	TITLE	PAGE No.
1.0	Pre-Approval	3
2.0	Objective	4
3.0	Scope	4
4.0	Responsibility	5
5.0	Equipment Details	6
6.0	Pre-Qualification Requirement	6
7.0	Tests & Checks	7
8.0	Checklist Of All Tests And Checks	14
9.0	Documents To Be Attached	15
10.0	Non Compliance	15
11.0	Deviation From Pre-Defined Specification, If Any	15
12.0	Change Control, If Any	15
13.0	Review Inclusive of Follow Up Action, If Any	16
14.0	Conclusion	16
15.0	Recommendations	16
16.0	Abbreviations	17
17.0	Post Approval	18



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

1.0 PRE – APPROVAL:

INITIATED BY:

DESIGNATION	NAME	SIGNATURE	DATE
OFFICER/EXECUTIVE (QUALITY ASSURANCE)			

REVIEWED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (PRODUCTION)			
HEAD (QUALITY CONTROL)			
HEAD (ENGINEERING)			

APPROVED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (QUALITY ASSURANCE)			



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

2.0 OBJECTIVE:

- To provide documented evidence that the Equipment is performing consistently, repeatedly and reproducibly within its established operating range and the results of all the test parameters meet the pre-defined acceptance criteria.
- To confirm the suitability of the Standard Operating Procedures for all routine activities associated with the system.

3.0 SCOPE:

- The scope of this report is limited for qualification of Vial Washing Machine installed in Washing and Sterilizing Room.
- This report provides all the relevant information of the performance qualification activity, In-process observations and analytical data of testing of collected samples.



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

4.0 RESPONSIBILITY:

The Validation Group, comprising of a representative from each of the following departments, shall be responsible for the overall compliance of this Protocol cum Report:

DEPARTMENTS	RESPONSIBILITIES
Quality Assurance	<ul style="list-style-type: none">• Preparation, Authorization, Approval and Compilation of the Performance Qualification.• Co-ordination with Quality Control, Production and Engineering to carryout Performance Qualification Activity.• Monitoring of Performance Qualification.
Production	<ul style="list-style-type: none">• Review of Protocol.• To co-ordinate and support Performance Qualification Activity.
Quality Control	<ul style="list-style-type: none">• Review of Protocol.• Analytical Support (Microbiological Testing/Analysis)
Engineering	<ul style="list-style-type: none">• Reviewing of qualification protocol for correctness, completeness and technical excellence• Responsible for trouble shooting (if occurred during execution).• Maintenance & preventive maintenance as per schedule.



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

5.0 EQUIPMENT DETAILS:

Equipment Name	Vial Washing Machine
Equipment
Manufacturer's Name	Ambica Pharma Machines Pvt. Ltd
Model	cGMP Model
Supplier's Name	Ambica Pharma Machines Pvt. Ltd
Location of Installation	Washing & Sterilization Room

6.0 PRE – QUALIFICATION REQUIREMENTS:

Verification for availability, completeness and approval status of all the required relevant documents shall be done and observations shall be recorded in the performance qualification report.

- Executed and approved Design Qualification document.
- Executed and approved Installation Qualification document.
- Executed and approved Operational Qualification document.
- SOP for Operation & Cleaning of Vial Washing Machine.
- SOP for Preventive Maintenance Vial Washing Machine.



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

7.0 TESTS AND CHECKS:

7.1 Verification of Documents:

Record the observations for documents in the below mentioned table.

S.No.	Document Name	Document/SOP No.	Completed (Yes/No)	Checked By (Engineering) Sign/Date	Verified By (QA) Sign/Date
1.	Executed and approved Design Qualification document				
2.	Executed and approved Installation Qualification document				
3.	Executed and approved Operational Qualification document				
4.	PQ Protocol approved				
5.	SOP for Operation & Cleaning of Vial Washing Machine				
6.	SOP for Preventive Maintenance Vial Washing Machine				

**Checked By
(Production)**

Sign/Date:

Verified By

(Quality Assurance)

Sign/Date:

Inference:

.....
.....
.....

**Reviewed By
(Manager QA)**

Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

7.2 Charcoal Test

Date of Test	
Vial Size	
Vial type	
Cycle	
Machine Speed	
Compressed Air Pressure	
WFI Pressure	
Re-circulated Water Pressure	
Purified Water Pressure	

Vial no.	Test	Observation			Analysed by Quality Control (Sign/Date)
		Initial	Middle	End	
1	VP				
2	VP				
3	VP				
4	VP				
5	VP				
6	VP				
7	VP				
8	VP				
9	VP				
10	VP				
11	VP				
12	VP				
13	VP				
14	VP				
15	VP				
16	VP				
17	VP				
18	VP				



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

Vial no.	Test	Observation			Analysed by Quality Control (Sign/Date)
		Initial	Middle	End	
19	VP				
20	VP				
	Average SVP for 20 Vials				

Acceptance criteria:

Visible particles (VP) : The vials should be free from visible particles.

Sub visible particles (SVP) : $\geq 10 \mu$: Not more than 6000

$\geq 25 \mu$: Not more than 600

**Checked By
(Production)**

Sign/Date:

Verified By

(Quality Assurance)

Sign/Date:

Inference:

.....
.....
.....
.....

Reviewed By

(Manager QA)

Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

7.3 Chloride Test

Date of Test	
Vial Size	
Vial type	
Cycle	
Machine Speed	
Compressed Air Pressure	
WFI Pressure	
Re-circulated Water Pressure	
Purified Water Pressure	

Vial no.	Observation			Analysed by Quality Control (Sign/Date)
	Initial	Middle	End	
1				
2				
3				
4				
5				
6				
7				
8				
9				
10				
11				



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

Vial no.	Observation			Analysed by Quality Control (Sign/Date)
	Initial	Middle	End	
12				
13				
14				
15				
16				
17				
18				
19				
20				

Acceptance criteria: All individual washed vials should not show turbidity or opalescence after adding reagents.

**Checked By
(Production)
Sign/Date:**

**Verified By
(Quality Assurance)
Sign/Date:**

Inference:

.....
.....
.....
.....
.....

**Reviewed By
(Manager QA)
Sign/Date:**



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

7.4 Glass Particle Test

Date of Test	
Vial Size	
Vial type	
Cycle	
Machine Speed	
Compressed Air Pressure	
WFI Pressure	
Re-circulated Water Pressure	
Purified Water Pressure	

Vial no.	Observation			Analysed by Quality Control (Sign/Date)
	Initial	Middle	End	
1				
2				
3				
4				
5				
6				
7				
8				
9				
10				



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

Vial no.	Observation			Analysed by Quality Control (Sign/Date)
	Initial	Middle	End	
11				
12				
13				
14				
15				
16				
17				
18				
19				
20				

Acceptance criteria: The vials should be free from the glass particles (Visual inspection).

**Checked By
(Production)**

Sign/Date:

Verified By

(Quality Assurance)

Sign/Date:

Inference:

.....

.....

.....

.....

.....

Reviewed By

(Manager QA)

Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

8.0 CHECKLIST OF ALL TESTS & CHECKS:

This checklist is provided to ensure that all tests or checks required for this protocol have been executed.

Tests or Checks	Executed (Yes/No)	Remarks
Verification of DQ, IQ & OQ & Other Documents		
Verification of Machine Performance		

**Checked By
(Production)
Sign/Date:**

**Verified By
(Quality Assurance)
Sign/Date.....**

Inference:

.....

.....

.....

.....

.....

**Reviewed By
(Manager QA)
Sign/Date:**



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

9.0 DOCUMENTS TO BE ATTACHED:

- Operation and Maintenance Manual.
- Copy of SOP's.
- Any Other Relevant Documents.

10.0 NON COMPLIANCE:

.....
.....
.....
.....
.....
.....

11.0 DEVIATION FROM PREDEFINED SPECIFICATION IF, ANY:

.....
.....
.....
.....
.....
.....
.....
.....
.....

12.0 CHANGE CONTROL, IF ANY:

.....
.....
.....
.....
.....
.....



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

13.0 REVIEW (INCLUSIVE OF FOLLOW UP ACTION, IF ANY):

.....
.....
.....
.....
.....
.....
.....

14.0 CONCLUSION:

.....
.....
.....
.....
.....
.....
.....
.....
.....

15.0 RECOMMENDATION:

.....
.....
.....
.....
.....
.....
.....
.....
.....



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

16.0 ABBREVIATIONS:

Sr.	:	Senior
Asst.	:	Assistant
No.	:	Number
WHO	:	World Health Organization
FDA	:	Food and Drug Administration
CFR	:	Code of Federal Regulations
cGMP	:	Current Good Manufacturing Practices
EU	:	European Union
QA	:	Quality Assurance
IQ	:	Installation Qualification
mm	:	Millimetre
Amp.	:	Ampere
DQ	:	Design Qualification
IQ	:	Installation Qualification
OQ	:	Operational Qualification
PQ	:	Performance Qualification
SOP	:	Standard Operating Procedure
Kg	:	Kilogram
No.	:	Number
Ltd.	:	Limited



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
VIAL WASHING MACHINE**

PROTOCOL No.:

17.0 POST APPROVAL:

INITIATED BY:

DESIGNATION	NAME	SIGNATURE	DATE
OFFICER/EXECUTIVE (QUALITY ASSURANCE)			

REVIEWED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (PRODUCTION)			
HEAD (QUALITY CONTROL)			
HEAD (ENGINEERING)			

APPROVED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (QUALITY ASSURANCE)			