



**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

**PERFORMANCE QUALIFICATION
REPORT
FOR
AUTOCARTONATOR**

EQUIPMENT ID. No.

LOCATION

PACKING HALL

DATE OF QUALIFICATION

SUPERSEDES PROTOCOL No.

Nil



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

REPORT CONTENTS

S.No.	TITLE	PAGE No.
1.0	Report Pre-Approval	3
2.0	Objective	4
3.0	Scope	4
4.0	Responsibility	5
5.0	Equipment Details	6
6.0	Pre-Qualification Requirement	6
7.0	Tests & Checks	7
8.0	Checklist Of All Tests And Checks	11
9.0	Documents To Be Attached	11
10.0	Non Compliance	12
11.0	Deviation From Pre-Defined Specification, If Any	12
12.0	Change Control, If Any	12
13.0	Review Inclusive of Follow Up Action, If Any	12
14.0	Conclusion	12
15.0	Recommendations	13
16.0	Abbreviations	13
17.0	Report Post-Approval	14



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

1.0 REPORT PRE – APPROVAL:

INITIATED BY:

DESIGNATION	NAME	SIGNATURE	DATE
OFFICER/EXECUTIVE (QUALITY ASSURANCE)			

REVIEWED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (PRODUCTION)			
HEAD (QUALITY CONTROL)			
HEAD (ENGINEERING)			

APPROVED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (QUALITY ASSURANCE)			



**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

2.0 OBJECTIVE:

- To provide documented evidence that the Equipment is performing consistently, repeatedly and reproducibly within its established operating range and the results of all the test parameters meet the pre-defined acceptance criteria.
- To confirm the suitability of the Standard Operating Procedures for all routine activities associated with the system.

3.0 SCOPE:

- The scope of this report is limited for qualification of Autocartonator installed in Packing Hall.
- This report provides all the relevant information of the performance qualification activity, In-process observations and analytical data of testing of collected samples.



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

4.0 RESPONSIBILITY:

The Validation Group, comprising of a representative from each of the following departments, shall be responsible for the execution of Performance Qualification Report.

DEPARTMENTS	RESPONSIBILITIES
Quality Assurance	<ul style="list-style-type: none">• Preparation, Pre-Approval and Compilation of the Performance Qualification Report.• Co-ordination with Quality Control, Production and Engineering to carryout Performance Qualification Activity.• Monitoring of Performance Qualification.• Post Approval of Performance Qualification Report after Execution.
Production	<ul style="list-style-type: none">• Review of Performance Qualification Report.• To co-ordinate and support Performance Qualification Activity.• Post Approval of Performance Qualification Report after Execution.
Quality Control	<ul style="list-style-type: none">• Analytical Support (Microbiological Testing/Analysis).
Engineering	<ul style="list-style-type: none">• Reviewing of qualification protocol for correctness, completeness and technical excellence• Responsible for trouble shooting (if occurred during execution).• Maintenance & preventive maintenance as per schedule.• Post Approval of Performance Qualification Report after Execution.



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

5.0 EQUIPMENT DETAILS:

Equipment Name	Autocartonator
Equipment	
Manufacturer's Name	Wimco ltd.
Supplier's Name	Wimco ltd.
Location of Installation	Packing Hall

6.0 PRE – QUALIFICATION REQUIREMENTS:

Verification for availability, completeness and approval status of all the required relevant documents shall be done and observations shall be recorded in the performance qualification report.

6.1 Verification of Documents:

Record the observations for documents in the below mentioned table.

S.No.	Document Name	Document/SOP No.	Completed (Yes/No)	Checked By (Engineering) Sign/Date	Verified By (QA) Sign/Date
1.	DQ Protocol approved				
2.	IQ Protocol approved				
3.	OQ Protocol approved				
4.	PQ Protocol approved				
5.	SOP for Operation & Cleaning of Autocartonator				
6.	SOP for Preventive Maintenance Autocartonator				

Checked By
Production
Sign/Date:

Verified By
Quality Assurance
Sign/Date:

Inference:

.....
.....
.....

Reviewed By
Manager QA
Sign/Date:



**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

7.0 TESTS AND CHECKS:

7.1 TRIAL NO.:

Date of test		Product Name		
Batch No.		Pack Size		
Speed		Low Speed (40 tubes/ min)		
Tube No.	Printing matter	Stereo Impression	Damage Carton	Scratches
1				
2				
3				
4				
5				
6				
7				
8				
9				
10				

RESULTS: The test results Complies / Not Complies as per Specification.

Checked By
Production
Sign/Date:

Verified By
Quality Assurance
Sign/Date:

Inference:

.....
.....
.....

Reviewed By
Manager QA
Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

7.2 TRIAL No.:

Date of test		Product Name	
Batch No.		Pack Size	

Speed	Optimum Speed (70 tubes/ min)
--------------	--------------------------------------

Tube No.	Printing matter	Stereo Impression	Damage Carton	Scratches
1				
2				
3				
4				
5				
6				
7				
8				
9				
10				

RESULTS: The test results Complies / Not Complies as per Specification.

Checked By
Production
Sign/Date:

Verified By
Quality Assurance
Sign/Date:

Inference:

.....
.....
.....

Reviewed By
Manager QA
Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

7.3 Trial No.:

Date of test		Product Name	
Batch No.		Pack Size	

Speed	High Speed (100 tubes/ min)
--------------	------------------------------------

Tube No.	Printing matter	Stereo Impression	Damage Carton	Scratches
1				
2				
3				
4				
5				
6				
7				
8				
9				
10				

RESULTS: The test results Complies / Not Complies as per Specification.

Checked By
Production
Sign/Date:

Verified By
Quality Assurance
Sign/Date:

Inference:

.....

Reviewed By
Manager QA
Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

7.4 Calculation of Rejection:

Trial No.	Speed (Tubes/min.)	No. of Tubes Checked	Passed Tubes	Rejected Tubes
1	40			
	70			
	100			
2	40			
	70			
	100			
3	40			
	70			
	100			
Total Tubes Taken`			Rejection %	
Rejection NMT 2.0 %				

Checked By
Production
Sign/Date:

Verified By
Quality Assurance
Sign/Date:

Inference:

.....
.....
.....

Reviewed By
Manager QA
Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

8.0 CHECKLIST OF ALL TESTS & CHECKS:

This checklist is provided to ensure that all tests or checks required for this protocol have been executed.

Tests or Checks	Executed (Yes/No)	Remarks
Verification of Documents		
Speed Synchronization		
Uniformity of Product		

Checked By
Production
Sign/Date:

Verified By
Quality Assurance
Sign/Date:.....

Inference:

.....

.....

.....

.....

.....

Reviewed By
Manager QA
Sign/Date:



PHARMA DEVILS

**PERFORMANCE QUALIFICATION REPORT
FOR
AUTOCARTONATOR**

PROTOCOL No.:

9.0 DOCUMENTS TO BE ATTACHED:

- Executed Raw Data.
- Any Other Relevant Documents.

10.0 NON COMPLIANCE:

.....
.....
.....
.....
.....

11.0 DEVIATION FROM PREDEFINED SPECIFICATION IF, ANY:

.....
.....
.....
.....
.....

12.0 CHANGE CONTROL, IF ANY:

.....
.....
.....
.....
.....

13.0 REVIEW (INCLUSIVE OF FOLLOW UP ACTION, IF ANY):

.....
.....
.....
.....
.....

14.0 CONCLUSION:

.....

.....

.....

.....

.....

.....

15.0 RECOMMENDATION:

.....

.....

.....


.....

.....

.....

16.0 ABBREVIATIONS:

- Asst. : Assistant
- cGMP : Current Good Manufacturing Practices
- CQA : Corporate Quality Assurance
- PQ : Performance Qualification
- Vol. : Volume
- i.e. : That is
- SS : Stainless steel
- Ltr. : Litre
- Nos. : Numbers.
- SOP : Standard Operating Procedure
- SS : Stain less Steel
- OFS : Autocartonator
- WHO : World Health Organization

 PHARMA DEVILS	PERFORMANCE QUALIFICATION REPORT FOR AUTOCARTONATOR	PROTOCOL No.:
--	--	----------------------

17.0 REPORT POST-APPROVAL:

INITIATED BY:

DESIGNATION	NAME	SIGNATURE	DATE
OFFICER/EXECUTIVE (QUALITY ASSURANCE)			

REVIEWED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (PRODUCTION)			
HEAD (QUALITY CONTROL)			
HEAD (ENGINEERING)			

APPROVED BY:

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (QUALITY ASSURANCE)			