

# PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE				
<b>Department:</b> Production (External Preparation)	SOP No.:			
Title: Operation and Cleaning of Air Jet Cleaning Machine	Effective Date:			
Supersedes: Nil	<b>Review Date:</b>			
Issue Date:	Page No.:			

## **1.0 OBJECTIVE:**

To lay down a procedure for Operation and Cleaning of Air jet cleaning machine.

### **2.0 SCOPE:**

This SOP is applicable for Operation and Cleaning of Air jet cleaning machine.

#### 3.0 **RESPONSIBILITY**:

Officer / Executive – Production

### 4.0 ACCOUNTABILITY:

Head – Production

#### 5.0 ABBREVIATIONS:

IPA	Iso Propyl Alcohol
Ltd.	Limited
MCB	Miniature Circuit Board
No.	Number
PO	Production Ointment
Pvt.	Private
QA	Quality Assurance
SOP	Standard Operating Procedure

### 6.0 **PROCEDURE**:

### 6.1 PRECAUTIONS AND CHECK POINTS:

- **6.1.1** Ensure that proper electric supply is provided to the machine.
- **6.1.2** Ensure the cleanliness of the machine, & surrounding area is free from any unwanted materials or balance materials of previous products.
- 6.1.3 Ensure that all removable panels of machine are at the specified place.
- 6.1.4 Check the pressure range of Compressed air,

### 6.2 MACHINE SETTING & OPERATION:

- **6.2.1** Before start the Air jet cleaning Machine set the required change parts.
- **6.2.2** Check the electrical main supply of Air jet cleaning Machine.
- **6.2.3** Start the main MCB (inside the panel).
- 6.2.4 Start main switch of Power Supply.



# PHARMA DEVILS

PRODUCTION DEPARTMENT

Department: Production (External Preparation)	SOP No.:
Title: Operation and Cleaning of Air Jet Cleaning Machine	<b>Effective Date:</b>
Supersedes: Nil	<b>Review Date:</b>
Issue Date:	Page No.:

# 6.3 **OPERATION:**

- **6.3.1** Start the De-Cartooning procedure in De-Cartooning Area as per SOP.
- 6.3.2 Transfer the bottle & Jar to inbuilt turntable of Air jet cleaning Machine.
- **6.3.3** Bottle transfer to Inverter-1 by Star wheel for cleaning.
- **6.3.4** After reaching the cleaning section air nozzles starts to flush the air inside towards the Bottle and simultaneously vacuum will suck the particles in Inverter-2, disturbed by the air.
- 6.3.5 Enter the operation details in Equipment Log in **Reference SOP.**

## 6.4 CLEANING:

- 6.4.1 Switch "OFF" the Machine.
- **6.4.2** Remove the side guard.
- **6.4.3** For stain mop the machine outer surface 70% IPA solution.
- 6.4.4 Make an entry of Cleaning Time in Equipment Log in **Reference SOP.**

### 7.0 ANNEXURES:

Not Applicable.

**ENCLOSURES:** SOP Training Record

## 8.0 **DISTRIBUTION:**

- Controlled Copy No. 01 Quality Assurance
- Controlled Copy No. 02 Production
- Master Copy
  Quality Assurance

### 9.0 **REFERENCES:**

Party Manual

### **10.0 REVISION HISTORY:**

# CHANGE HISTORY LOG

Revision No.	Change control No.	Details of Changes	Reason for Change	Effective Date	Updated By