PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE

Title: Procedure for Charging of Vials, Dropper and Cap into the Hopper

SOP No.:		Department:	Production
SOP No.:		Effective Date:	
Revision No.:	00	Revision Date:	
Supersede Revision No.:	Nil	Page No.:	1 of 2

1.0 **OBJECTIVE**:

To lay down a procedure for charging of vials, dropper and cap into the hopper.

2.0 SCOPE:

This SOP is applicable for charging of vials, dropper and cap into the hopper of vials filling, capping and screwing machine of Three Piece Line.

3.0 **RESPONSIBILITY:**

Officer / Executive Production

4.0 **ACCOUNTABILITY:**

Head Production

5.0 ABBREVIATIONS:

IPA Iso Propyl Alcohol LAF Laminar Air Flow QA Quality Assurance

SOP Standard Operating Procedure

Pvt. Private Ltd. Limited

6.0 PROCEDURE:

6.1 INSTRUCTIONS:

- **6.1.1** Sanitized the hands with filtered 70% IPA Solution before touching the bag containing vials, dropper and cap.
- **6.1.2** Take the bags of vial, dropper and cap placed under LAF in buffer room and transfer it into filling room by using the mobile LAF.
- 6.1.3 Take out the bags of vial, dropper and cap from mobile LAF and put it on the table of the bench type LAF.
- 6.1.4 Take out the bags of vial, dropper and cap from bench type LAF and transfer it into the Filling LAF and open the bag aseptically with the help of knife blade.
- **6.1.5** Then transfer the vial, dropper and cap in to respective hopper.
- **6.1.6** Place the loose bag of vials, dropper and cap under Bench LAF aseptically.

7.0 ANNEXURES:

Not Applicable

ENCLOSURES: SOP Training Record



PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE

Title: Procedure for Charging of Vials, Dropper and Cap into the Hopper

SOP No.:		Department:	Production
SOP No.:	Effective Date:		
Revision No.:	00	Revision Date:	
Supersede Revision No.:	Nil	Page No.:	2 of 2

8.0 DISTRIBUTION:

• Controlled Copy No.01 Quality Assurance

• Controlled Copy No.02 Production

• Master Copy Quality Assurance

9.0 REFERENCES:

Not Applicable.

10.0 REVISION HISTORY:

CHANGE HISTORY LOG

Revision	Change	Details of Changes	Reason for Change	Effective	Updated
No.	Control No.			Date	$\mathbf{B}\mathbf{y}$