

PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE

Title: Loading of Bottles in Autoclave Trolley

SOP No.:		Department:	Production
SOF No.:		Effective Date:	
Revision No.:	00	Revision Date:	
Supersede Revision No.:	Nil	Page No.:	1 of 2

1.0 OBJECTIVE:

To lay down a Procedure for Loading of Filled / Sealed Bottles in Autoclave Trolley.

2.0 SCOPE:

This SOP is applicable for storage of loading of filled /sealed bottles in autoclave trolley in Production Area.

3.0 RESPONSIBILITY:

Officer / Executive Production

4.0 ACCOUNTABILITY:

Head Production

5.0 ABBRIVATION:

BPCR Batch Production & Control Record

Ltd Limited


LVP Large Volume Parenteral

No. Number PVT Private

SOP Standard Operating procedure

6.0 PROCEDURE:

- **6.1** Loading of filled /sealed bottles in autoclave trolley:
 - **6.1.1** Collect the filled and sealed containers of LVP products coming out of the filling area in cleaned perforated trays of a trolley.
 - **6.1.2** Arrange the 330 bottles (15X 22) in a tray.
 - **6.1.3** Arrange as per given diagram:

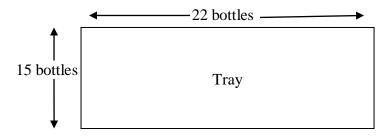

- **6.1.4** Load 12 trays in a trolley (330 X 12) -3960 bottles in a trolley.
- **6.1.5** Record the quantity in respective log and BPCR.

6.1.6 PRECAUTIONS:



PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE

Title: Loading of Bottles in Autoclave Trolley

SOP No.:		Department:	Production
SOF No.:		Effective Date:	
Revision No.:	00	Revision Date:	
Supersede Revision No.:	Nil	Page No.:	2 of 2

- **6.1.6.1** Before starting loading of bottles of a new product / batch, check that no any bottles of previous product/ batch are lying on a tray/ trolly.
- **6.1.6.2** Affix the chemical indicator strip/3m tape to each trolley before loading in to autoclave & affix after unloading the batch in respective print out.

7.0 ANNEXURES:

Not Applicable

ENCLOSURES: SOP Training Record

8.0 DISTRIBUTION:

Controlled Copy No.01
Quality Assurance

• Controlled Copy No.02 Production

• Master Copy Quality Assurance

9.0 REFERENCES:

Not Applicable.

10.0 REVISION HISTORY:

CHANGE HISTORY LOG

Revision No.	Change control No.	Details of Changes	Reason for Change	Effective Date	Updated By