

PROTOCOL No.:

# PERFORMANCE QUALIFICATION

### **REPORT**

### **FOR**

### **COMMINUTING MILL**

EQUIPMENT ID No.	
LOCATION	Granulation
DATE OF QUALIFICATION	
SUPERSEDES REPORT No.	NIL



PROTOCOL No.:

### **CONTENTS**

S.No.	TITLE	PAGE No.
1.0	Report Pre-Approval	3
2.0	Objective	4
3.0	Scope	4
4.0	Responsibility	5
5.0	<b>Equipment Details</b>	6
6.0	Pre-Qualification Requirements	6
7.0	Tests & Checks	7-8
8.0	Checklist of all Tests & Checks	9
9.0	Documents Attached	9
10.0	Non Compliance	9
11.0	Deviation From Pre-Defined Specification, If Any	10
12.0	Change Control, If Any	10
13.0	Review (Inclusive of follow up action, If Any)	10
14.0	Conclusion	10
15.0	Recommendation	10
16.0	Abbreviations	11
17.0	Report Post Approval	12



P	R	O'	T	()	$\mathbf{C}$	$\Omega$	Γ,	N	O.	•

\L:

PREPARED BY:

DESIGNATION	NAME	SIGNATURE	DATE
OFFICER/EXECUTIVE (QUALITY ASSURANCE)			

### **REVIEWED BY:**

DESIGNATION	NAME	SIGNATURE	DATE
OPERATING MANAGER (QUALITY ASSURANCE)			
HEAD (ENGINEERING)			

### **APPROVED BY:**

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (PRODUCTION)			



PROTOCOL No.:

#### 2.0 OBJECTIE:

• To provide documented evidence that the Equipment is performing consistently, repeatedly and reproducibly within its established operating range and the results of all test parameters meet the pre-defined acceptance criteria

#### 3.0 SCOPE:

- This Report covers all the aspects of Performance Qualification for the **Comminuting Mill** (Make: .....) installed in **Granulation**.
- This Report will define the methods and documentation used to qualify the Comminuting Mill for PQ.

.



PROTOCOL No.:

### 4.0 **RESPONSIBILITY:**

The Validation Group, comprising of a representative from each of the following departments, shall be responsible for the overall compliance of this Report.

DEPARTMENTS	RESPONSIBILITIES
	Preparation, Authorization, review and Compilation of the
	Performance Qualification.
<b>Quality Assurance</b>	<ul> <li>Co-ordination with Quality Control, Production and Engineering to</li> </ul>
	carryout Performance Qualification Activity.
	Monitoring of Performance Qualification.
Production	Approval of Report.
Troduction	• To co-ordinate and support Performance Qualification Activity.
	Reviewing of qualification Report for correctness, completeness and
Engineering	technical excellence
Engineering	• Responsible for trouble shooting (if occurred during execution).
	Maintenance & preventive maintenance as per schedule.



PROTOCOL No.	PRO	TO	COL	No.
--------------	-----	----	-----	-----

5.0 EQUIPMENT DETAIL	PMENT DETAIL	S
----------------------	--------------	---

<b>Equipment Name</b>	Comminuting Mill
Equipment ID.	
Model	
Sr. No.	
Manufacturer's Name	Cadmach Machinery Pvt. Ltd.
Supplier's Name	Cadmach Machinery Pvt. Ltd.
<b>Location of Installation</b>	Granulation

### **6.0 PRE – QUALIFICATION REQUIREMENTS:**

### **6.1 Verification of Documents:**

Record the observations for documents in the below mentioned table.

S.No.	Document Name	Document /SOP No.	Completed (Yes/No)	Checked By (Engineering) Sign/Date	Verified By (Quality Assurance) Sign/Date
1.	PQ Protocol approved				
2.	SOP for Operation & Cleaning of Comminuting Mill				
3.	SOP for Preventive  Maintenance Comminuting  Mill				

Inference:	
	Reviewed By
	(Manager QA)
	(Manager QA) Sign & Date:



PRO	TO	COL	No.:
$\mathbf{I}$	$\mathbf{U}$	COL	110

#### 7.0 TESTS AND CHECKS:

#### 7.1 TEST PRODUCT BATCH INFORMATION:

S.No.	Product Name	Batch No.	Batch Size	Mfg. Date	Expiry				
Drug Products									
Compil (QA) Sign/D	led By ate:								
Inferen	nce:								
		•••••	•••••		•••••				
•••••		•••••••	••••••	•••••	•••••				
••••••		••••••	•••••	•••••	•••••				
			$(\mathbf{N}$	eviewed By Ianager QA) gn/Date:					



**Acceptance Criteria** 

### PERFORMANCE QUALIFICATION REPORT FOR COMMINUTING MILL

### 7.2 REPORT OF PERFORMANCE EVALUATION USING DRUG PRODUCTS:

#### **In-Process Observations:**

Product Name/	Screen specified	Sieve specified	% Granules	% Granules	Black	Integrity of Screen & sieve	
Batch No.	as per BMR	as per BMR	Passed	Retained	Particles	Before	After

% Retain Granules	NMT 5%
Black Particles	Should be absent
Integrity of Screen & Sieve	Should be Integrated
Checked By	Verified By
(Production)	(Quality Assurance)
Sign/Date:	Sign/Date:
Inference:	
	Reviewed By
	(Manager QA)
	Sign/Date:



P	R	O'	T	()	$\mathbf{C}$	$\Omega$	Γ,	N	O.	•

### 8.0 CHECKLIST OF ALL TESTS & CHECKS:

This checklist is provided to ensure that all tests or checks required for this protocol have been executed.

Executed

	Tests or Checks	(Yes/No)	Remarks
Verific	ation of Performance using Drug Products.		
Checke	d By		Verified By
(Produc			(Quality Assurance)
Sign/Da	te:		Sign/Date
Inferen	ce:		
• • • • • • • • • • • • • • • • • • • •			
• • • • • • • • • • • • • • • • • • • •			
			Reviewed By (Manager QA)
			Sign/Date:
9.0 D	OOCUMENTS TO BE ATTACHED:		
•	Any other relevant document.		
0.0 N	ON COMPLIANCE:		



n	-	$\sim$		$\sim$	C	$\sim$	r	<b>T</b> 4		
ν									Λ	•
	77	v	1	v	v	C,	┙.	Τ.	v.	•

11.0	DEVIATION FROM PRE-DEFINED SPECIFICATION, IF ANY:
12.0	CHANGE CONTROL, IF ANY:
13.0	REVIEW (INCLUSIVE OF FOLLOW UP ACTION, IF ANY)
14.0	CONCLUSION
15.0	RECOMMENDATION



PROTOCOL No.:

#### **16.0 ABBREVIATIONS:**

CDL : Comminuting Mill

IQ : Installation Qualification

Kg : Kilogram
Ltd : Limited

MOC : Material of Construction

NLT : Not Less Than

PQ : Performance Qualification

Pvt : Private

QA : Quality Assurance

SOP : Standard Operating Procedure



P	R	$\cap$	T(	$\mathbf{C}$	$\mathbf{O}$	Vo.	•

### 17.0 REPORT POST – APPROVAL:

#### PREPARED BY:

DESIGNATION	NAME	SIGNATURE	DATE
OFFICER/EXECUTIVE (QUALITY ASSURANCE)			

### **REVIEWED BY:**

DESIGNATION	NAME	SIGNATURE	DATE
OPERATING MANAGER (QUALITY ASSURANCE)			
HEAD (ENGINEERING)			

### **APPROVED BY:**

DESIGNATION	NAME	SIGNATURE	DATE
HEAD (PRODUCTION)			