



PHARMA DEVILS

MICROBIOLOGY DEPARTMENT

STANDARD OPERATING PROCEDURE

Title: Cleaning and Operation of Vacuum Pump

SOP No.:		Department:	Microbiology
		Effective Date:	
Revision No.:	00	Revision Date:	
Supersede Revision No.:	Nil	Page No.:	1 of 2

1.0 OBJECTIVE

To lay down procedure for cleaning and operation of Vacuum pump.

2.0 SCOPE

This SOP is applicable for all the Vacuum pumps, Make - Millipore used in Quality control Department.

3.0 RESPONSIBILITY

Prepared by - Executive Quality Control

Checked by - Assistant Manager Microbiology / QC

Approved by - Head QA, QC

4.0 PROCEDURE

4.1 Operation

- 4.1.1 Connect the vacuum pump plug the main supply.
- 4.1.2 Connect the vacuum pump to the one side of the filtrate collecting bottle through silicon tubing.
- 4.1.3 Connect the other side of the filtrate-collecting bottle to the micro fill assembly through silicon tubing.
- 4.1.4 Switch on the main supply.
- 4.1.5 Start the vacuum pump by pressing the button provided on the vacuum pump.
- 4.1.6 Operate the procedure till filtration is complete.
- 4.1.7 After completion of filtration switch off the vacuum pump by pressing the button.
- 4.1.8 Switch off the main supply.
- 4.1.9 Remove the silicon tubing from the filtrate-collecting bottle and vacuum pump.

4.2 Cleaning

- 4.2.1 Insure that the power supply to the vacuum pump is switched off and main cord is removed from the supply.
- 4.2.2 Wipe the vacuum pump with a clean cloth wetted with disinfectant solution everyday.

5.0 SAFETY & PRECAUTIONS

Before cleaning or any maintenance remove the power supply cord of the vacuum pump from the main supply.

6.0 REVISION HISTORY



PHARMA DEVILS

MICROBIOLOGY DEPARTMENT

STANDARD OPERATING PROCEDURE

Title: Cleaning and Operation of Vacuum Pump

SOP No.:		Department:	Microbiology
		Effective Date:	
Revision No.:	00	Revision Date:	
Supersede Revision No.:	Nil	Page No.:	2 of 2

Revision No.	Reason for Revision	Superseded from & date
00	First Issue	-----

7.0 REFERENCES

Not applicable.

8.0 ABBREVIATIONS

SOP : Standard Operating Procedure

9.0 ANNEXURES

Not applicable