



PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE

Department: Production

SOP No.:

Title: Cleaning of Leak Test Apparatus

Effective Date:

Supersedes: Nil

Review Date:

Issue Date:

Page No.:

Purpose: To lay down operating procedure for the cleaning of Leak Test Apparatus.

Scope: This procedure is applicable for

Responsibility: Production Chemist
To supervise the activity

Pre-operating Check: Check no packing material or residue of previous product should be there.

Frequency: For different product each time after use or at the end of the shift for the same product.

Equipment used: Leak test apparatus

Cleaning procedure:

1. Discard the water from the apparatus chamber
2. Clean the plate, lid etc with DM.WATER.
3. Mop the above with clean cloth then that no residue of product there.
4. Mop the desicator and m/c and also clean the place where the Equipment is kept and it's surrounding area.