



PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE

Department: Production

SOP No.:

Title: Cleaning of Stirrer

Effective Date:

Supersedes: Nil

Review Date:

Issue Date:

Page No.:

Purpose: To lay down operating procedure for the Cleaning of Stirrer.

Scope: This procedure is applicable for

Responsibility: Production Chemist
To supervise the activity

Pre-operating Check: Check no material or residue should be there.

Frequency: For different product each time after use or at the end of the shift for the same product.

Equipment used: Stirrer

Cleaning procedure:

1. Unscrew the propeller and rinse the rode and propeller with portable water and finally with DM Water.
2. Clean the propeller and rod with brush to remove any residue of the previous product and then wash with portable water and finally with DM. water.
3. Clean the stirrer with dry and clean cloth.