



PHARMA DEVILS

PRODUCTION DEPARTMENT

STANDARD OPERATING PROCEDURE

Department: Production	SOP No.:
Title: SOP for Operation of Dust Extractor	Effective Date:
Supersedes: Nil	Review Date:
Issue Date:	Page No.:

Purpose: To lay down operating procedure for the cleaning of Dust Extractor.

Scope: This procedure is applicable to

Responsibility: Production Chemist
To supervise the activity

Pre-operating Check: Ensure that the line clearance is done.

Equipment used: Dust extractor

Operating procedure:

1. Tie the finger in the dust extractor.
2. Fit the s.s. Container (storage device) in dust extractor.
3. Fit the vacuum pipe of dust extractor to the dust collection clamp.
4. Switch on the dust extractor.
5. At the end of shift or after completion of tablet compression removed the dust powder collected in the s.s.container (storage device) and destroyed the powder.