



# PHARMA DEVILS

PRODUCTION DEPARTMENT

## STANDARD OPERATING PROCEDURE

<b>Department:</b> Production	<b>SOP No.:</b>
<b>Title:</b> SOP for Operation of Tablet Deduster	<b>Effective Date:</b>
<b>Supersedes:</b> Nil	<b>Review Date:</b>
<b>Issue Date:</b>	<b>Page No.:</b>

**Purpose:** To lay down operating procedure for the operation of tablet deduster

**Scope:** This procedure is applicable for .....

**Responsibility:** Production Chemist  
To supervise the activity

**Pre-operating Check:** Check no material or residue should be there.

**Equipment used:** Tablet deduster

### Wet cleaning procedure:

1. Check the cleaning of deduster and surrounding area.
2. Assemble the all parts of tablet deduster.
3. Tie the dust outlet pipe with polythylene bag.
4. Set the tablet deduster below the tablet outlet channel of compression.
5. Plug the tablet deduster into suitable power supply point and switch on.
6. After completion of tablet compression removed and destroyed dust powder collected in polythylene bag.