

PHARMA DEVILS

QUALITY ASSURANCE DEPARTMENT

System ID:	Document Number:	
Effective Date: Version Number: 00		
eam Quality Test		Test Run: _
		he chamber is o
 Steam quality test kit Sampling Bottles Tube SOP for steam quality testing 		
Equipment should be normal operation	nal mode	
More t	han Test Description	
Check the quality of steam for non conde	nsable gases as per the SOP	
Check the quality of steam for superheat	gases as per the SOP	
Check the quality of steam for dryness va	lue as per the SOP	
Record the observations & data as per Ap	ppendix 3.1.1.	
Acceptance	criteria	Acceptance criteria fulfilled? (Y/N
The fraction of non-condensable gases sh	ould not exceed 3.5%.	
The superheat measured in the expansion	tube should not exceed 25°C	
	-	
Comm	ent	Deviation Ref.
1	eam Quality Test Objective if this test is to demonstrate to requisite physical, chemical & microbiologous Steam quality test kit Sampling Bottles Tube SOP for steam quality testing Equipment should be normal operations More to the Check the quality of steam for non conder the Check the quality of steam for dryness vans Record the observations & data as per Appearance The fraction of non-condensable gases should not be less that be within ±3°C of the value measured during the steam of the condensable gases of the value measured during the condensable gases of the value measured during the condensable gases of the condensable gases	te: Version Number: 00 eam Quality Test Objective if this test is to demonstrate that the steam being introduced into t requisite physical, chemical & microbiological quality. • Steam quality test kit • Sampling Bottles • Tube