



VALIDATION PROTOCOL FOR MIXING OF INGREDIENTS IN PLANETARY MIXER

1. Objective:

To validate the process of Dry Mixing of Raw materials so as to establish that the blend uniformity of active ingredient is achieved at standard dry mixing time and the blend uniformity results are within the specified limit.

2. Scope:

Applicable to all blending/lubrication operation in PLM.

3. Principle:

Mixing of raw materials due to rotation of Agitator in Planetary Mixer.

4. Site of the Study:

Hormone Department.

5. Responsibility:

Production :

Quality Control :

Engineering :

Quality Assurance:

(Individuals to be named in the report)

6. Description of the Equipment to be used:

PLANETARY MIXER: λ

CODE No. : Details to be recorded in the report.

CAPACITY : Details to be recorded in the report.

RPM (STIRRER) : Details to be recorded in the report.

λ Date of Equipment Qualification to be recorded in the Report.

Sampling thief and Bulk sampler Code No. to be recorded in the validation report.

7. BMR and Standard Operating Procedure (SOP) to be followed:

- i) SOP for operating Planetary Mixer: SOP number to be recorded in the report.
- ii) SOP for sampling with sampling thief: SOP number to be recorded in the report.
- iii) Batch Manufacturing Record: Formulation Code No. and Manufacturing code No.: To be recorded in the validation report

8. Control:

8.1 Requirement:

- i. Active Pharmaceutical Ingredient to be used should meet the requirement of Specification (Name of the active ingredient, Quantity required, Pharmacopoeia grade and the



VALIDATION PROTOCOL FOR MIXING OF INGREDIENTS IN PLANETARY MIXER

test involved for raw material testing with result and Analytical Reference number should be recorded in Validation report).

- ii. Validated Analytical Methods for estimation of active ingredient (Reference Analytical Validation protocol number / Reference Specification No. to be recorded in the report).
- iii. Analytical Reference number for validation Technical Information Sheet (T. I. Sheet).

8.2 Calibration:

Calibration details of PLM (Calibration details to be recorded in the report).

8.3 Training:

Availability of Training Record of Personnel involved in the Validation Exercise should be recorded in the Validation Report

8.4 Precaution:

Safety aspects while operation of equipment and process must be ensured.

9. Validation Procedure:

- 9.1 Load the sifted ingredients in the Planetary Mixer as per the sequence given in the Batch Manufacturing Record (BMR).
- 9.2 Operate the Planetary Mixer as per SOP.
- 9.3 Sample one to three unit doses in triplicate from 10 positions at various depths as shown in sampling plan on Page 6, with the help of a sampling thief, at standard mixing time specified in the BMR. Send these samples to Quality Control for Content Analysis.
- 9.4 The results should be reported as Individual test results and calculate % RSD.
- 9.5 Check the appearance of mix after completion of mixing operation and record the same in report.

10. Acceptance criteria:

The Optimal time decided should conform to the following:

- 10.1 Uniform distribution of the mix, when checked visually.
- 10.2 Content of Active Ingredient upon testing as per Quality Control Specification should be within the specified limit.
- 10.3 The Relative Standard Deviation (RSD) of content of Active Ingredient sampled at the optimal time should be less than 5%.

11. Non Compliances:



VALIDATION PROTOCOL FOR MIXING OF INGREDIENTS IN PLANETARY MIXER

Details of deviations (including justification of acceptance) done to successfully carry out the validation exercise and any OOS results obtained should be checked. (Attach the details in the Validation report).

12. Type of Validation:

Concurrent Validation/Re-validation

13. Frequency:

- 1) Concurrent validation : Three consecutive successful validation exercises.
- 2) Re-validation (Periodic) : One validation exercise within five years.
- 3) Revalidation (after major change) : Three consecutive successful validation exercises.

14. Results/Observations:

Record the observations during the study and results obtained from Quality Control Department in the Validation Report.

15. Summary of findings of experiment (inference):

Summarize the findings of the Validation Study to draw an inference.

16. Recommendation (Including requirements of any additional Documentation):

Record the recommendations based on the interpretation of the results of the Validation Report.

17. Team approval:

The individuals who have performed the Validation Study, supervised the validation, completed the records, performed the testing of the product should approve the validation report.

18. Review (inclusive of follow up action, if any):

The Validation Report should be reviewed by Unit Quality Assurance and Unit Head. The report should include any follow up action if required.

19. Approved by:

Unit Quality Assurance and Unit Head should finally approve the validation Report.

20. Attachments:

Annexure (if any) attached to the Validation Report should be recorded.

20. Abbreviations:

VP : Validation Protocol



PHARMA DEVILS

QUALITY ASSURANCE DEPARTMENT

VALIDATION PROTOCOL FOR MIXING OF INGREDIENTS IN PLANETARY MIXER


Lt. : Litres
No. : Number
A.R.No. : Analytical Reference Number
RSD : Relative Standard Deviation
mg : Milligram
OOS : Out of Specification
% : Percent
SOP : Standard Operating Procedure
BMR : Batch Manufacturing Record
RPM : Revolution Per Minutes




VALIDATION PROTOCOL FOR MIXING OF INGREDIENTS IN PLANETARY MIXER

SAMPLING PLAN FOR BLENDING/LUBRICATION

TOP VIEW OF PLANETARY MIXER (PLM)

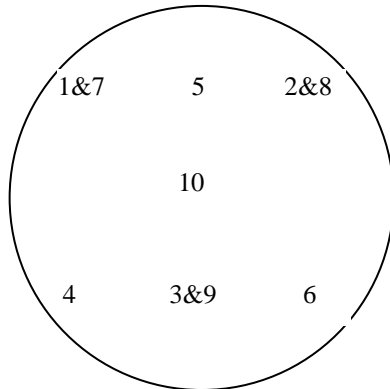

SIDE VIEW OF PLANETARY MIXER

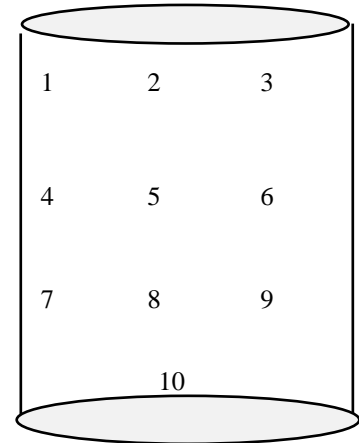

SAMPLING POSITIONS:

POSITION 1, 2 & 3 : Top Layer of the bed.

POSITION 4, 5 & 6 : 0.5 feet from top layer of the bed

POSITION 7, 8 & 9 : 1 feet from top layer of the bed

POSITION 10 : Bottom centre where the tip touches the PLM



PHARMA DEVILS

QUALITY ASSURANCE DEPARTMENT

VALIDATION PROTOCOL FOR MIXING OF INGREDIENTS IN PLANETARY MIXER